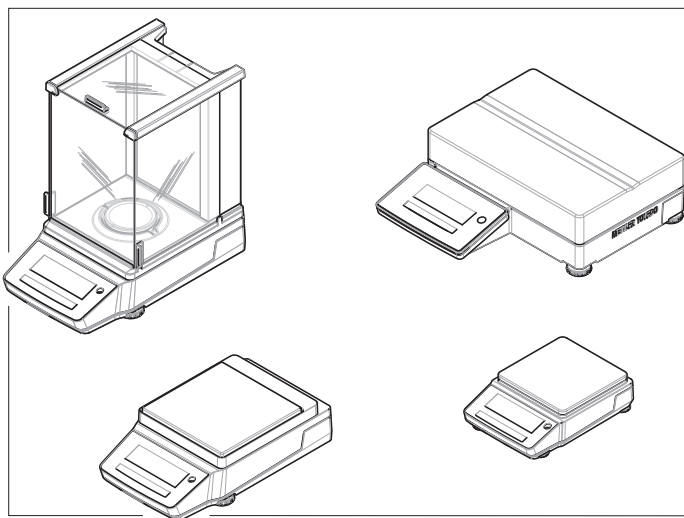
English	User Manual Analytical and Precision Balances MA
Deutsch	Benutzerhandbuch Analysen- und Präzisionswaagen MA
Español	Manual de usuario Balanzas analíticas y de precisión MA
Français	Guide de l'utilisateur Balances d'analyse et de précision fine MA
简体中文	用户手册 分析与精密天平 MA
日本語	ユーザマニュアル 分析天びん・上皿天びん MA
Čeština	Návod k použití Analytické a přesné váhy MA
Dansk	Brugervejledning Analyse- og præcisionsvægte MA
Hrvatski	Korisnički priručnik Analitičke i precizne vage MA
Magyar	Felhasználói útmutató Analitikai és precíziós mérlegek MA

METTLER TOLEDO

► www.mt.com/MA-RM

en For full information, always refer to the Reference Manual (RM).

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User Manual **Analytical and Precision Balances**

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Magyar

Table of Contents

1	Introduction	3
1.1	Document purpose	3
1.2	Further documents and information	3
1.3	Acronyms and abbreviations	3
2	Safety Information	4
2.1	Definition of signal words and warning symbols	4
2.2	Product-specific safety notes	4
3	Design and Function	5
3.1	Overview analytical balances	6
3.2	Overview precision balances, small	6
3.3	Overview precision balances, large	7
3.4	Overview precision balances, compact	7
3.5	Overview terminal	8
3.6	Overview interface connections	9
3.7	Components description	9
3.7.1	Draft shield	9
3.7.2	Weighing pan	9
3.7.3	Drip tray	9
3.7.4	Door handle	10
3.7.5	Leveling feet	10
3.7.6	Terminal	10
3.7.7	QuickLock for doors and front panel	10
3.7.8	Release button for back panel	11
3.8	User interface	11
3.8.1	Main sections at a glance	11
3.8.2	Main weighing screen	12
4	Installation and Putting into Operation	14
4.1	Selecting the location	14
4.2	Unpacking the balance	15
4.3	Installation	15
4.3.1	Balances with draft shield	16
4.3.2	Balances without draft shield	18
4.3.3	Balances, large	18
4.3.4	Balances, compact	18
4.4	Putting into operation	19
4.4.1	Connecting the balance	19
4.4.2	Switching on the balance	19
4.4.3	Leveling the balance	19
4.4.3.1	Leveling large balances	20
4.4.3.2	Leveling compact balances	20
4.4.3.3	Leveling all other balances	21
4.4.4	Performing an internal adjustment	21
4.4.5	Entering / Exiting standby mode	22
4.4.6	Entering / Exiting hibernation mode	22
4.4.7	Switching off the balance	22
4.5	Performing a simple weighing	22
4.5.1	Opening and closing draft shield doors	22
4.5.2	Zeroing the balance	22
4.5.3	Taring the balance	23
4.5.4	Performing a weighing	23

4.6	Using batteries (compact balance).....	23
4.6.1	Inserting or replacing batteries.....	23
4.7	Transporting, packing, and storing.....	24
4.7.1	Transporting the balance over short distances.....	24
4.7.2	Transporting the balance over long distances.....	24
4.7.3	Packing and storing.....	24
5	Maintenance	25
5.1	Maintenance tasks.....	25
5.2	Cleaning.....	25
5.2.1	Disassembling for cleaning.....	25
5.2.1.1	Balances with draft shield.....	26
5.2.1.2	Balances without draft shield.....	28
5.2.1.3	Balances, large.....	28
5.2.1.4	Balances, compact.....	29
5.2.2	Cleaning the balance.....	29
5.2.3	Putting into operation after cleaning.....	30
5.3	Service.....	30
6	Technical Data	30
6.1	General data.....	30
7	Disposal	31
8	Compliance Information	31

1 Introduction

Thank you for choosing a METTLER TOLEDO balance. The balance combines high performance with ease of use.

EULA

The software in this product is licensed under the METTLER TOLEDO End User License Agreement (EULA) for Software.

When using this product you agree to the terms of the EULA.

▶ www.mt.com/EULA

1.1 Document purpose

This User Manual provides brief instructions about the first steps to take with the instrument. This ensures safe and efficient handling. Personnel must have carefully read and understood this manual before performing any task.

1.2 Further documents and information

This document is available in other languages online.

▶ www.mt.com/MA-UM

Product page:

▶ www.mt.com/MA-balances

Instructions for cleaning a balance, "8 Steps to a Clean Balance":

▶ www.mt.com/lab-cleaning-guide

Search for software:

▶ www.mt.com/labweighing-software-download

Search for documents:

▶ www.mt.com/library

For further questions, please contact your authorized METTLER TOLEDO dealer or service representative.

▶ www.mt.com/contact

1.3 Acronyms and abbreviations

Original term	Explanation
AC	Alternating Current
ASTM	American Society for Testing and Materials
DC	Direct Current
EMC	Electromagnetic Compatibility
FCC	Federal Communications Commission
GWP	Good Weighing Practice
HID	Human Interaction Device
ID	Identification
IP	Ingress Protection
LAN	Local Area Network
LED	Light-Emitting Diode
LPS	Limited Power Source
MAC	Media Access Control

MT-SICS	METTLER TOLEDO Standard Interface Command Set
NA	Not Applicable
OIML	Organisation Internationale de Métrologie Légale (International Organization of Legal Metrology)
RAM	Random Access Memory
RM	Reference Manual
SOP	Standard Operating Procedure
UM	User Manual
USB	Universal Serial Bus
USP	United States Pharmacopeia

2 Safety Information

Two documents named "User Manual" and "Reference Manual" are available for this instrument.

- The User Manual is available online in various languages.
- A printed version of the User Manual is delivered with the instrument.
- The Reference Manual is available online. This manual contains a full description of the instrument and its use.
- Keep both documents for future reference.
- Include both documents if you transfer the instrument to other parties.

Only use the instrument according to the User Manual and the Reference Manual. If you do not use the instrument according to these documents or if the instrument is modified, the safety of the instrument may be impaired and Mettler-Toledo GmbH assumes no liability.

2.1 Definition of signal words and warning symbols

Safety notes contain important information on safety issues. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results. Safety notes are marked with the following signal words and warning symbols:

Signal words

DANGER	A hazardous situation with high risk, resulting in death or severe injury if not avoided.
WARNING	A hazardous situation with medium risk, possibly resulting in death or severe injury if not avoided.
CAUTION	A hazardous situation with low risk, resulting in minor or moderate injury if not avoided.
NOTICE	A hazardous situation with low risk, resulting in damage to the instrument, other material damage, malfunctions and erroneous results, or loss of data.

Warning symbols

General hazard

Notice

2.2 Product-specific safety notes

Intended use

This instrument is designed to be used by trained staff. The instrument is intended for weighing purposes.

Any other type of use and operation beyond the limits of use stated by Mettler-Toledo GmbH without consent from Mettler-Toledo GmbH is considered as not intended.

Responsibilities of the instrument owner

The instrument owner is the person holding the legal title to the instrument and who uses the instrument or authorizes any person to use it, or the person who is deemed by law to be the operator of the instrument. The instrument owner is responsible for the safety of all users of the instrument and third parties.

Mettler-Toledo GmbH assumes that the instrument owner trains users to safely use the instrument in their workplace and deal with potential hazards. Mettler-Toledo GmbH assumes that the instrument owner provides the necessary protective gear.

Safety notes

WARNING

Death or serious injury due to electric shock

Contact with parts that carry a live current can lead to death or injury.

- 1 Only use the METTLER TOLEDO power cable and AC/DC adapter designed for your instrument.
- 2 Connect the power cable to a grounded power outlet.
- 3 Keep all electrical cables and connections away from liquids and moisture.
- 4 Check the cables and the power plug for damage and replace them if damaged.

NOTICE

Damage to the instrument or malfunction due to the use of unsuitable parts

- Only use parts from METTLER TOLEDO that are intended to be used with your instrument.

A list of spare parts and accessories can be found in the Reference Manual.

3 Design and Function

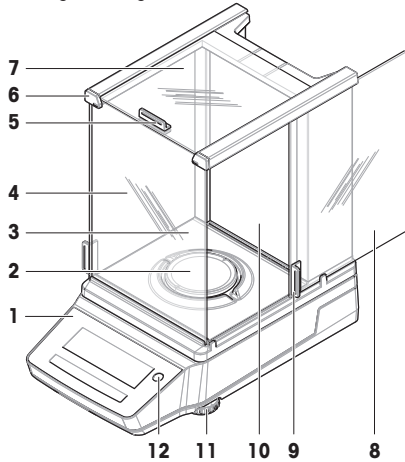
For further information, consult the Reference Manual (RM).

► www.mt.com/MA-RM

3.1 Overview analytical balances


0.01 mg / 0.1 mg


1	Terminal	7	Top door, draft shield
2	Weighing pan	8	Side door, draft shield (right/left)
3	Drip tray	9	Handle, side door
4	Front panel, draft shield	10	Back panel, draft shield
5	Handle, top door	11	Leveling feet
6	QuickLock, front panel	12	Level indicator

3.2 Overview precision balances, small

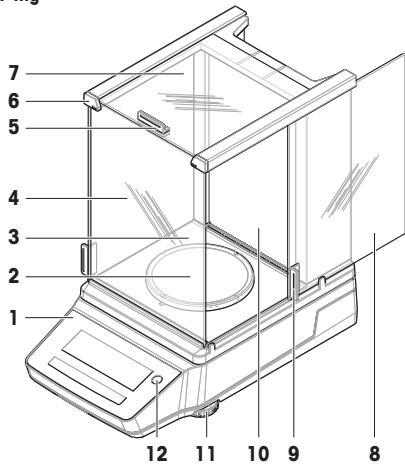1 mg


10 mg / 100 mg

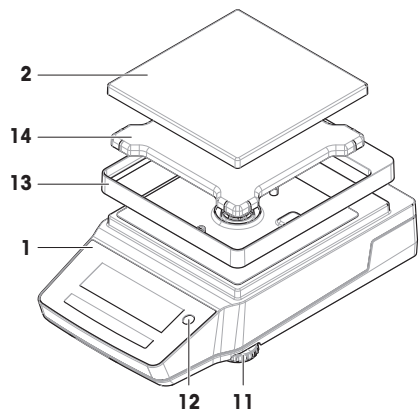1	Terminal	8	Side door, draft shield (right/left)
2	Weighing pan	9	Handle, side door
3	Drip tray	10	Back panel, draft shield
4	Front panel, draft shield	11	Leveling feet
5	Handle, top door	12	Level indicator
6	QuickLock, front panel	13	Draft-protection element
7	Top door, draft shield	14	Weighing pan support

3.3 Overview precision balances, large


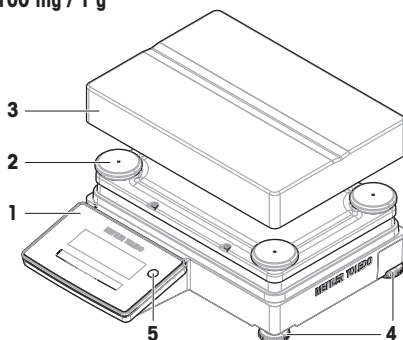
100 mg / 1 g

1	Terminal	4	Leveling feet
2	Weighing pan support cap	5	Level indicator
3	Weighing pan		

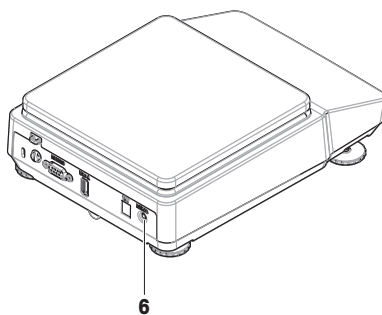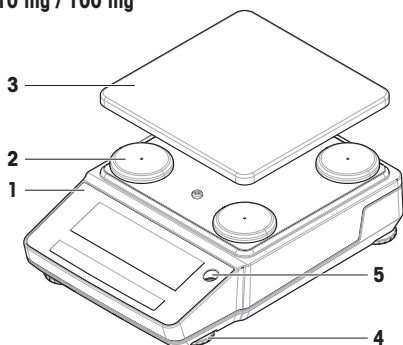
3.4 Overview precision balances, compact


10 mg / 100 mg

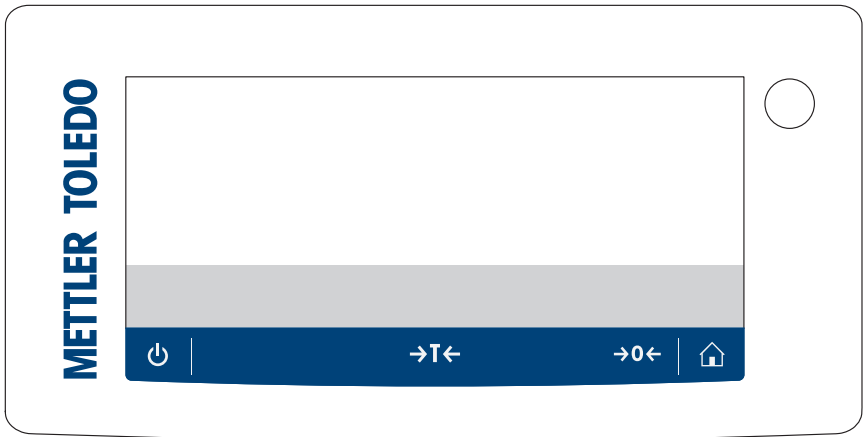
1	Terminal	4	Leveling feet
2	Weighing pan support cap	5	Level indicator
3	Weighing pan	6	Wake-up switch (battery mode)

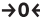3.5 Overview terminal

	Name	Description
	Standby	<p>By tapping , the balance is not completely switched off but goes into standby mode. To switch the balance completely off, it must be unplugged from the power supply.</p> <p>Note Do not disconnect the balance from the power supply unless the balance is not used for an extended period of time. After switching on the instrument, it must warm up before giving accurate results.</p> <p>Note Compact balances in battery mode only: By tapping , the balance is completely switched off. There is no standby mode.</p>
	Tare	<p>Tares the balance.</p> <p>This function is used when the weighing process involves containers. After taring the balance, the screen shows <i>Net</i> which indicates that all displayed weights are net.</p>
	Zero	<p>Zeroes the balance.</p> <p>The balance must always be zeroed before starting the weighing process. After zeroing, the balance sets a new zero point.</p>
	Home	<p>To return from any settings menu level to the main weighing screen of the currently set or last used application.</p>

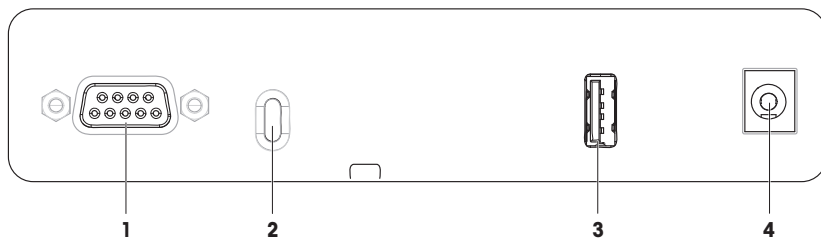3.6 Overview interface connections


1	RS232C serial interface	3	USB-A port
2	Slot for anti-theft cable	4	Socket for AC/DC adapter

3.7 Components description


3.7.1 Draft shield

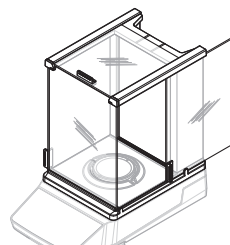The draft shield protects the weighing area against environmental impacts like drafts or moisture. The side doors and the top door can be opened manually.


3.7.2 Weighing pan

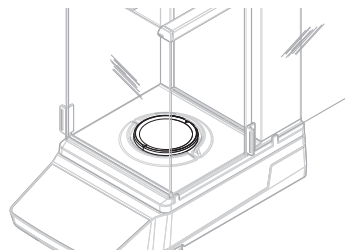The weighing pan is the load receptor that serves to accommodate the weighing item.


3.7.3 Drip tray

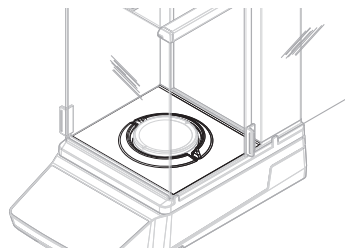The drip tray is positioned below the weighing pan. The primary purpose of the drip tray is to ensure quick cleaning of the balance.

3.7.4 Door handle

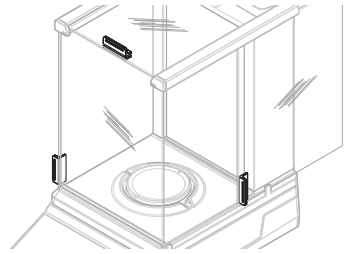The door handles are mounted on the draft shield doors. The handles are used to manually open the side doors and the top door of the draft shield.


3.7.5 Leveling feet

The balance stands on height-adjustable feet. These feet are used to level the balance.


Note

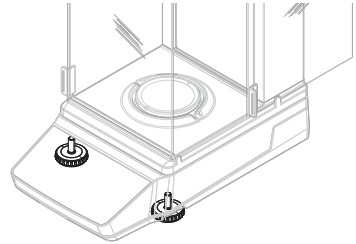Large balances and compact balances have four leveling feet.


3.7.6 Terminal

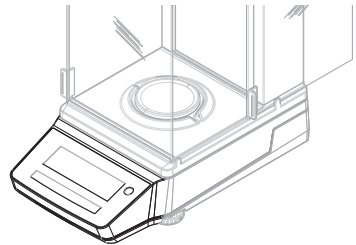The terminal is integrated into the balance and has an LCD display. The terminal and the weighing platform are protected by a replaceable cover.


3.7.7 QuickLock for doors and front panel

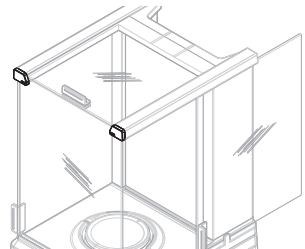Depending on the position, the QuickLock is used to lock/unlock the top door, the side doors, and the front panel of the draft shield.

3.7.8 Release button for back panel

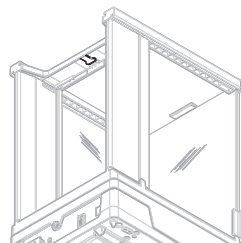The release button is used to lock/unlock the back panel of the draft shield.

3.8 User interface

3.8.1 Main sections at a glance

The main weighing screen (1) displays weighing results and provides feedback on the status of the balance. The operation keys (2) are the main navigation element where all the menus and settings can be accessed and selected.

	Name	Description
1	Main weighing screen	Displays the weighing results and provides information on the status of the balance.
2	Operation keys	Provides access to balance menus and functions.

Operation keys


	Name	Description
	Settings	Accesses the settings menu.
	Applications	Accesses the applications menu.
	Previous / Up / Increase	<ul style="list-style-type: none"> Scrolls up in a list of topics. Changes between unit 1, unit 2, and the application unit (if available). Increases numbers. Adds a captured weight within an application.
	Next / Down / Decrease	<ul style="list-style-type: none"> Scrolls down in a list of topics. Changes between unit 1, unit 2, and the application unit (if available). Decreases numbers.
	Cancel	Cancels a task or leaves the menu without saving.

	Name	Description
✓	Accept	<ul style="list-style-type: none"> • Accesses the selected menu. • Starts the selected application. • Confirms the entry.
	Adjustment	Executes the predefined adjustment procedure.
	Print / Transfer	<ul style="list-style-type: none"> • Prints the displayed value. • Prints the settings of the active user menu. • Transfers data to the predefined target location.

See also

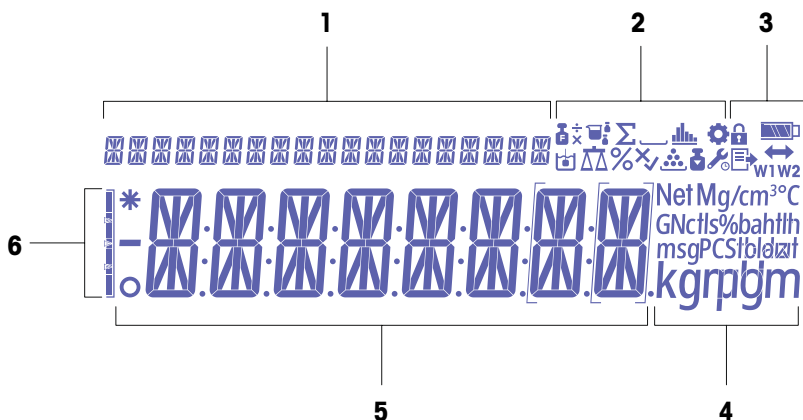Main weighing screen ▶ Page 12

3.8.2 Main weighing screen

1	Coach text	4	Unit icons
2	Application icons	5	Weight value / menu label / parameter label / parameter value
3	Status icons	6	Weighing aid

Application icons

	Name	Description
	Application "Division factor weighing"	The application "Division factor weighing" divides a predefined factor by the weight value (in grams): result = factor / weight The result is rounded to a predefined number of decimal places.
	Application "Multiplication factor weighing"	The application "Multiplication factor weighing" multiplies the weight value (in grams) by a predefined factor: result = factor × weight The result is rounded to a predefined number of decimal places.

	Name	Description
	Application "Formulation"	The application "Formulation" (Net total) has several purposes. <ul style="list-style-type: none"> • Weigh in (add and store) up to 999 individual component weights and display the total. If a printer is connected, the component weights are printed individually and as a total. • Tare/pre-tare and store up to 999 container weights and display the total. If a printer is connected, the tare weights are printed individually and as a total. • Fill up the sum of all component net weight values by adding a further component to a higher value.
	Application "Totaling"	The application "Totaling" allows you to weigh in different samples to add their weight values and to totalize them. 1 to 999 samples are possible.
	Application "Statistics"	The application "Statistics" allows you to generate statistics of a series of weighing values. 1 to 999 values are possible.
	Application "Dynamic weighing"	The application "Dynamic weighing" allows you to determine the weights of unstable samples or to determine weights under unstable ambient conditions. The balance calculates the weight as the average of a number of weighing operations over a defined time.
	Application Density	The application Density allows you to determine the density of solid bodies and liquids. Determination of the density uses "Archimedes' principle" according to which a body immersed in a fluid undergoes an apparent loss in weight which is equal to the weight of the fluid it displaces.
	Application Weighing	The application Weighing is used for simple weighing tasks, or to perform a measurement series.
	Application Check weighing	The application Check weighing allows you to check the deviation of a sample weight within a tolerance limit to a reference target weight.
	Application Counting	The application Counting allows you to determine the number of pieces put on the weighing pan.

Status icons

	Name	Description
	Menu settings activated	Menu settings are activated and can be adapted.
	Menu settings blocked	Menu settings are locked and cannot be adapted.
	Internal adjustment started	The internal adjustment of the balance has started.
	Service reminder	The next service is due.
	Publishing	Data is being published.
	Status of battery charge	Compact balances only: Indicates the level of the battery charge. <ul style="list-style-type: none"> • : Battery full • : Battery 3/4 full • : Battery 1/2 full • : Battery 1/4 full • : Battery discharged
	Connectivity	Data connection to another device.

	Name	Description
W1	Weighing range 1	Weighing range 1 is active.
W2	Weighing range 2	Weighing range 2 is active.
Net	Net weight values	"Net" indicates that all weight values displayed are net values.
G	Gross weight values	Gross weight values are displayed.
M	Stored value (Memory)	A stored value from the memory of the balance is displayed.

Label Icons

	Name	Description
*	Calculated value	Indicates that the displayed value is calculated.
o	Unstable value	Indicates the displayed value is unstable, meaning that it changes over time.
{ }	Uncertified digits	Brackets indicate uncertified digits (approved balance models only).

4 Installation and Putting into Operation

4.1 Selecting the location

A balance is a sensitive precision instrument. The location where it is placed will have a profound effect on the accuracy of the weighing results.

Requirements of the location

Place indoors on stable table

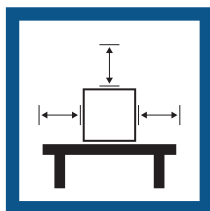
Ensure sufficient spacing

Level the instrument

Provide adequate lighting table


Avoid direct sunlight



Avoid vibrations

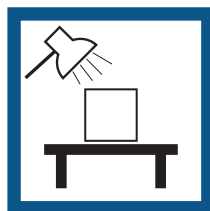Avoid strong drafts

Avoid temperature fluctuations

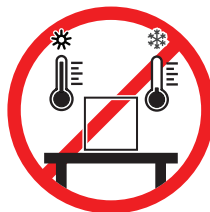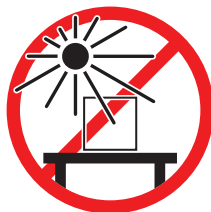Take into account the environmental conditions. See "Technical Data".

Sufficient spacing for balances: > 15 cm all around the instrument

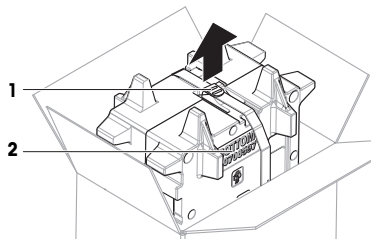
4.2 Unpacking the balance

Check the package, the packaging elements and the delivered components for damages. If any components are damaged, please contact your METTLER TOLEDO service representative.


Note

Depending on the balance model, the packaging elements and the components may look different.


- 1 Open the box and lift the package out using the lifting strap (1).
- 2 Place the package on a level surface with the inscription **BOTTOM (2)** facing downwards.

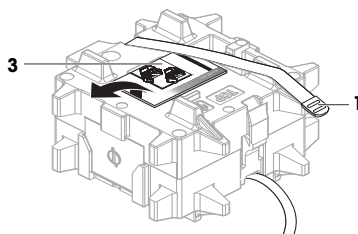
- 3 Open the lifting strap (1) and remove the User Manual (3).

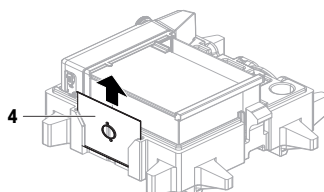- 4 Remove the upper part of the package and carefully unpack the drip tray (4).

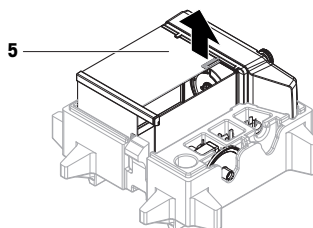- 5 Carefully unpack the balance (5) and all other items.
- 6 Remove the protective bag.
- 7 Keep the protective cover installed on the weighing platform and on the terminal.
- 8 Store all parts of the packaging in a safe place for future use.
 - ➔ The balance is ready for installation.

4.3 Installation

Note

Depending on the balance model, the components may look different.

4.3.1 Balances with draft shield

CAUTION

Injury due to sharp objects or broken glass

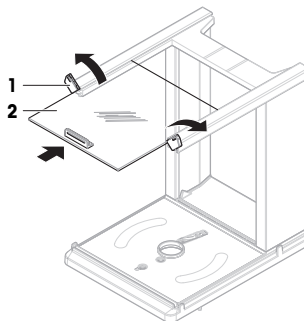
Instrument components, e.g., glass, can break and lead to injuries.

- Always proceed with focus and care.


Note

Skip steps 1 – 5 when unpacking the balance for the first time, or if the draft shield is already assembled and mounted on the weighing platform.


- 1 Assemble the draft shield: Turn the QuickLock (1, right, left) and slide in the top door (2).

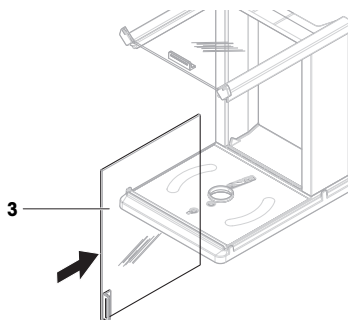
- 2 Slide in the side door (3) (right, left).

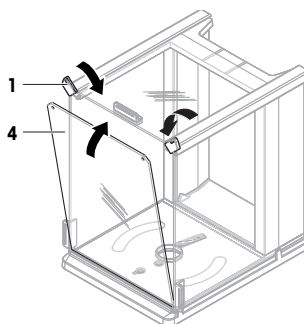
- 3 Attach the front panel (4), then turn the QuickLock (1, right, left) to hold the panel in place.

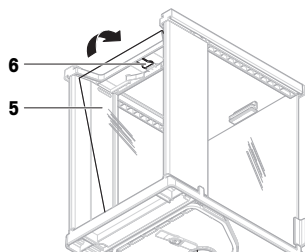
- 4 Attach the back panel (5). Make sure the release button (6) snaps in.

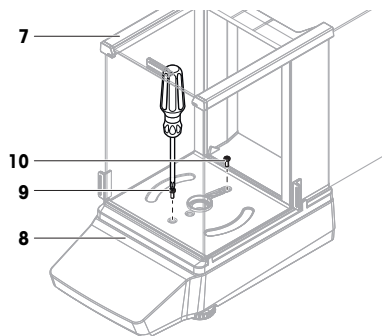
- 5 Secure the draft shield (7) to the weighing platform (8) by fixing the front screw (9) and the rear screw (10) with a Phillips screwdriver.

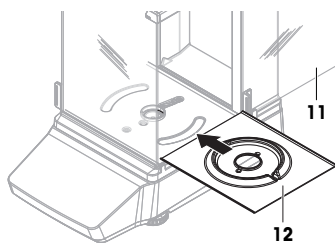- 6 Fully open the side door (11)
7 Insert the drip tray (12).

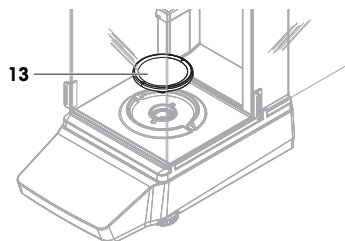
- 8 Install weighing pan (13).
➔ The balance is ready for use.


4.3.2 Balances without draft shield

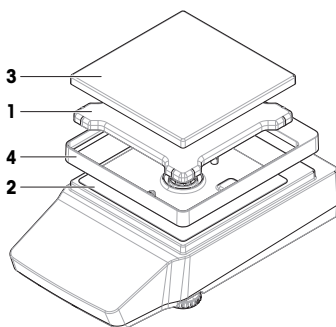- 1 Place the weighing pan support (1) on top of the weighing platform (2).
- Note**
To protect your balance, keep the protective cover installed on the weighing platform (2).
- 2 Place the weighing pan (3) on top of the weighing pan support (1).
- 3 Place the draft-protection element (4) on top of the weighing platform (2).
➔ The balance is ready for use.


4.3.3 Balances, large

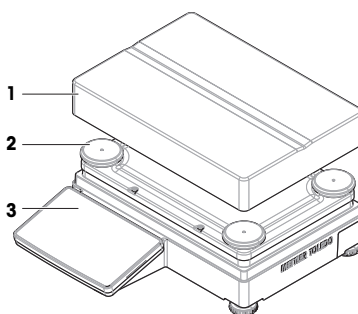- Place the weighing pan (1) on top of the support caps (2).
- Note**
To protect your balance, keep the protective cover installed on the terminal (3).
- ➔ The balance is ready for use.

4.3.4 Balances, compact

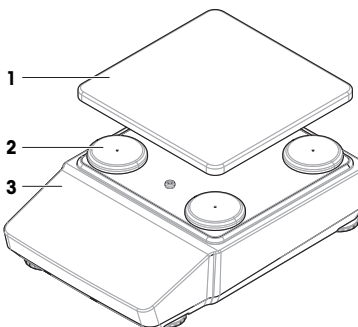- Place the weighing pan (1) on top of the support caps (2).
- Note**
To protect your balance, keep the protective cover installed on the weighing platform (3).
- ➔ The balance is ready for use.

4.4 Putting into operation

4.4.1 Connecting the balance

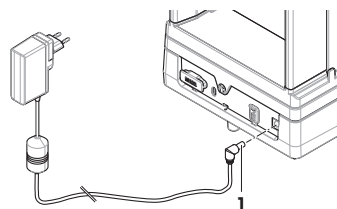
WARNING

Death or serious injury due to electric shock

Contact with parts that carry a live current can lead to death or injury.

- 1 Only use the METTLER TOLEDO power cable and AC/DC adapter designed for your instrument.
- 2 Connect the power cable to a grounded power outlet.
- 3 Keep all electrical cables and connections away from liquids and moisture.
- 4 Check the cables and the power plug for damage and replace them if damaged.

- 1 Install the cables in such a way that they cannot be damaged or interfere with operation.
- 2 Insert the plug of the AC/DC adapter (1) into the power socket of the instrument.
- 3 Insert the plug of the power cable into a grounded power outlet that is easily accessible.
 - ➔ The balance automatically switches on.

Note

Do not connect the instrument to a power outlet controlled by a switch. After switching on the instrument, it must warm up before giving accurate results.

See also

🔗 General data ▶ Page 30

4.4.2 Switching on the balance

When connected to the power supply, the balance automatically switches on.

Acclimatization and warm up

Before the balance gives reliable results, it must:

- acclimatize to the room temperature
- warm up by being connected to the power supply

The acclimatization time and warm-up time for balances are available in "General data".

Note

When the balance is exiting standby, it is ready immediately.

See also

🔗 General data ▶ Page 30

🔗 Entering / Exiting standby mode ▶ Page 22

🔗 Entering / Exiting hibernation mode ▶ Page 22

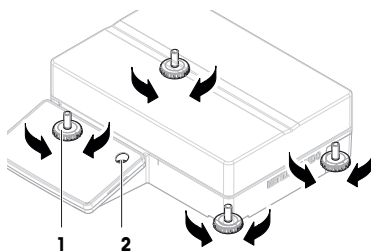
🔗 Switching off the balance ▶ Page 22

4.4.3 Leveling the balance

Exact horizontal and stable positioning are essential for repeatable and accurate weighing results.

4.4.3.1 Leveling large balances

- 1 Screw in all leveling feet fully.
- 2 Screw out all leveling feet two to three turns except for the rear left leveling foot.
- 3 Turn the leveling feet (1) until the air bubble is in the center of the level indicator (2). Proceed as shown in the following example.
- 4 Screw out the rear left leveling foot until it makes contact with the table surface.
 - ➔ The balance is leveled and supported by all four leveling feet.

Example

Air bubble at 12 o'clock:

- Turn both front feet clockwise.
- Turn rear right foot counterclockwise.

Air bubble at 3 o'clock:

- Turn left front foot clockwise, turn right front foot counterclockwise.
- Turn rear right foot counterclockwise.

Air bubble at 6 o'clock:

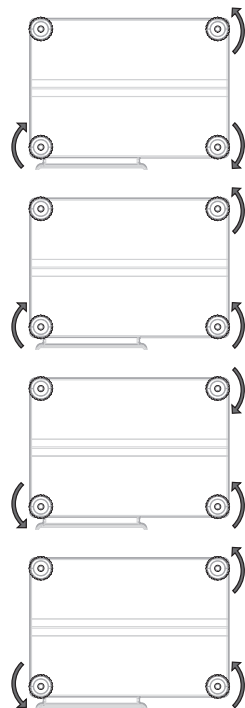
- Turn both front feet counterclockwise.
- Turn rear right foot clockwise.

Air bubble at 9 o'clock:


- Turn both front feet counterclockwise.
- Turn rear right foot counterclockwise.

4.4.3.2 Leveling compact balances

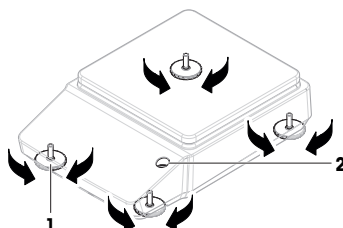- 1 Screw in both rear leveling feet fully.
- 2 Screw out both front leveling feet two to three turns.
- 3 Turn the front leveling feet (1) until the air bubble is in the center of the level indicator (2). Proceed as shown in the following example.
- 4 Screw out both rear leveling feet until they make contact with the table surface.
 - ➔ The balance is leveled and supported by all four leveling feet.

Example

Air bubble at 12 o'clock:

Turn both feet clockwise.

Air bubble at 3 o'clock:

Turn left foot clockwise, turn right foot counterclockwise.

Air bubble at 6 o'clock:

Turn both feet counterclockwise.


Air bubble at 9 o'clock:

Turn left foot counterclockwise, turn right foot clockwise.

**4.4.3.3 Leveling all other balances**

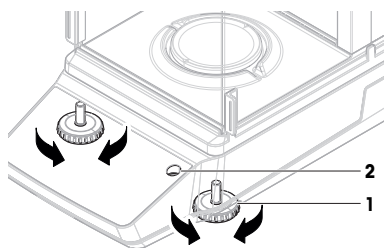
- Turn the leveling feet (1) until the air bubble is in the center of the level indicator (2).

**Example**

Air bubble at 12 o'clock:

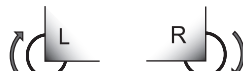
Turn both feet clockwise.

Air bubble at 3 o'clock:

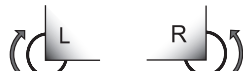Turn left foot clockwise, turn right foot counterclockwise.

Air bubble at 6 o'clock:

Turn both feet counterclockwise.

Air bubble at 9 o'clock:

Turn left foot counterclockwise, turn right foot clockwise.

**4.4.4 Performing an internal adjustment**

- 1 On the main weighing screen, press .
 - ➔ ADJ.INT appears.
- 2 Press .
 - ➔ The adjustment is executed.
 - ➔ The adjustment results appear.
- 3 Press .
 - ➔ The balance is ready.

4.4.5 Entering / Exiting standby mode

- 1 To enter standby mode, press  for less than 2 s.
 - ➔ The brightness of the display is reduced, information on the display is visible. The balance is still switched on.
- 2 To exit standby mode, press .
Alternatively, tap the weighing pan, or place a weight on the weighing pan.
 - ➔ The display is turned on.

4.4.6 Entering / Exiting hibernation mode

 **Note**

Only relevant for compact balances in battery mode.

- 1 To enter hibernation mode, press  for less than 2 s.
 - ➔ The display is turned off. The balance is in hibernation mode.
- 2 To exit hibernation mode, press the wake-up switch on the rear side of the balance.
 - ➔ The display is turned on.

4.4.7 Switching off the balance

To completely switch off the balance, it must be disconnected from the power supply. By pressing , the balance goes only into standby mode.

 **Note**

When the balance has been completely switched off for some time, it must warm up before it can be used.

See also

-  Switching on the balance ▶ Page 19
-  Entering / Exiting standby mode ▶ Page 22
-  Entering / Exiting hibernation mode ▶ Page 22

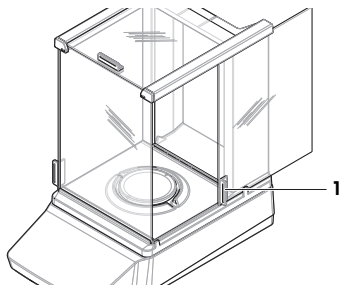
4.5 Performing a simple weighing

 **Note**

A balance with draft shield is used to explain the procedure. For balances without a draft shield, skip the instructions steps concerning the draft shield.

4.5.1 Opening and closing draft shield doors

- Open the door manually with the door handle (1).


4.5.2 Zeroing the balance

- 1 Open the draft shield.
- 2 Clear the weighing pan.
- 3 Close the draft shield.
- 4 Press  to zero the balance.
 - ➔ The balance is zeroed.

4.5.3 Taring the balance

If a sample vessel is used, the balance must be tared.

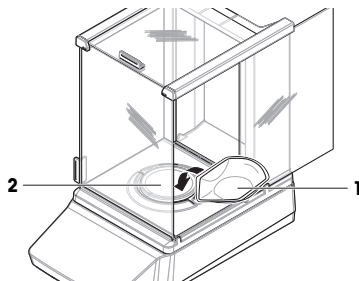- The balance is zeroed.
- 1 Place the sample vessel (1) on the weighing pan (2).
 - 2 Press →T← to tare the balance.
 - ⇒ The balance is tared. The icon **Net** appears.

4.5.4 Performing a weighing

- 1 Open the draft shield.
- 2 Place the weighing object into the sample vessel.
- 3 Close the draft shield.
- 4 Wait until the instability detector **O** disappears.
 - ⇒ The result is displayed.
- 5 Optional, if a printer is connected: Press  to print the weighing result.

4.6 Using batteries (compact balance)

The balance can also be operated with batteries. Under normal operating conditions, the balance runs independently of the AC power supply for about 8 to 15 hours (using alkaline batteries).

Immediately after an interruption of the AC power supply, the balance automatically switches to battery operation. This can happen, for example, by pulling the power plug, or in the event of a power failure. After the AC power supply is restored, the balance automatically switches back to AC operation.

The balance uses eight standard AA batteries (LR6). Alkaline batteries are preferred.

Rechargeable batteries can be used. Charging batteries inside the balance is **not** possible.

When the balance is operated with batteries, the battery symbol in the display lights up. The number of segments that are shown in the battery symbol indicate the charge level. When the batteries are almost completely discharged, the battery symbol flashes.

See also

 Main weighing screen ▶ Page 12

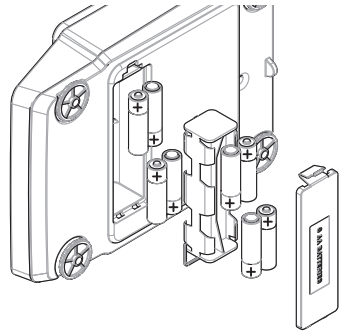
4.6.1 Inserting or replacing batteries

- Read and follow all warnings and instructions supplied by the battery manufacturer.
- Do not mix different types or brands of batteries. Performance of batteries varies depending on the manufacturer.
- Remove the batteries from the balance if the balance is not used for a long period of time.
- Batteries must be disposed of properly, according to local regulations.

Proceed as follows:

- The balance is switched off.
 - The weighing pan is removed.
- 1 Turn the balance carefully on its side.

- 2 Open and remove the battery-chamber cover.
- 3 Insert / replace the batteries with the correct polarity as shown in the battery holder.
- 4 Insert and close the battery-chamber cover.
- 5 Turn the balance carefully to its normal position.
- 6 Reinstall all components in the reverse order.


4.7 Transporting, packing, and storing

4.7.1 Transporting the balance over short distances

- 1 Disconnect the AC/DC adapter and unplug all interface cables.
- 2 Hold the balance with both hands and carry it in horizontal position to the target location. Consider the requirements of the location.

If you want to put the balance into operation, proceed as follows:

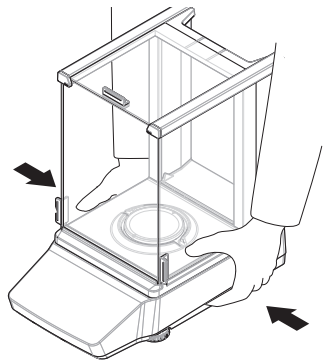- 1 Connect in reverse order.
- 2 Give the balance sufficient time to warm up.
- 3 Level the balance.
- 4 Perform an internal adjustment.

See also

- 🔗 Selecting the location ▶ Page 14
- 🔗 Switching on the balance ▶ Page 19
- 🔗 Leveling the balance ▶ Page 19
- 🔗 Performing an internal adjustment ▶ Page 21

4.7.2 Transporting the balance over long distances

METTLER TOLEDO recommends using the original packaging for transportation or shipment of the balance or balance components over long distances. The elements of the original packaging are developed specifically for the balance and its components and ensure maximum protection during transportation.

See also

- 🔗 Unpacking the balance ▶ Page 15

4.7.3 Packing and storing

Packing the balance

Store all parts of packaging in a safe place. The elements of the original packaging are developed specifically for the balance and its components, and ensures maximum protection during transportation and storage.

Storing the balance

Only store the balance under the following conditions:

- Indoor and in the original packaging

- According to the environmental conditions, see "Technical Data"

Note

When storing for longer than 6 months, the rechargeable battery may become empty (only date and time get lost).

See also

 Technical Data ▶ Page 30

5 Maintenance

To guarantee the functionality of the balance and the accuracy of the weighing results, a number of maintenance actions must be performed by the user.

For further information, consult the Reference Manual (RM).

www.mt.com/MA-RM

5.1 Maintenance tasks

Maintenance action	Recommended interval	Remarks
Performing an internal adjustment	<ul style="list-style-type: none"> • Daily • After cleaning • After leveling • After changing the location 	see "Performing an internal adjustment"
Performing routine tests (eccentricity test, repeatability test, sensitivity test). METTLER TOLEDO recommends to at least perform a sensitivity test.	<ul style="list-style-type: none"> • After cleaning • After assembling the balance • After a software update • Depending on your internal regulations (SOP) 	see "Tests" in the Reference Manual
Cleaning	<ul style="list-style-type: none"> • After every use • Depending on the degree of pollution • Depending on your internal regulations (SOP) 	see "Cleaning"
Updating the software	<ul style="list-style-type: none"> • Depending on your internal regulations (SOP). • After a new software release. 	see "Software update" in the Reference Manual

See also

 Performing an internal adjustment ▶ Page 21

 Cleaning ▶ Page 25

5.2 Cleaning

5.2.1 Disassembling for cleaning

Note

Depending on the balance model, the components may look different.

Note

In most cases, it is not necessary to remove the protective cover to clean the balance.

5.2.1.1 Balances with draft shield

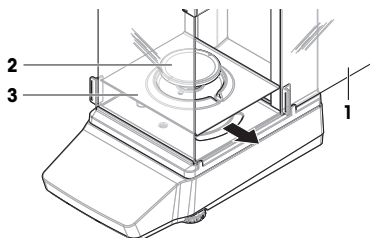
CAUTION

Injury due to sharp objects or broken glass


Instrument components, e.g., glass, can break and lead to injuries.

- Always proceed with focus and care.


- 1 Fully open the side door (1).
- 2 Remove the weighing pan (2) and the drip tray (3).

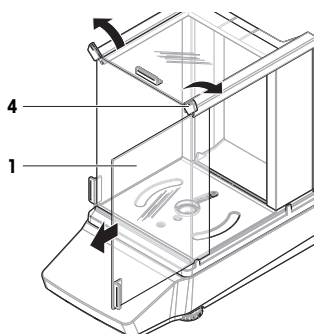
- 3 Turn the quick lock (4) and pull the side door (1) towards the front to remove it (right, left).

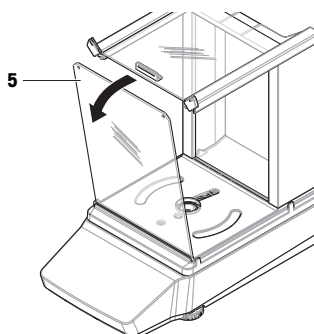- 4 Tilt the front panel (5) towards the front and lift it upwards to remove it.

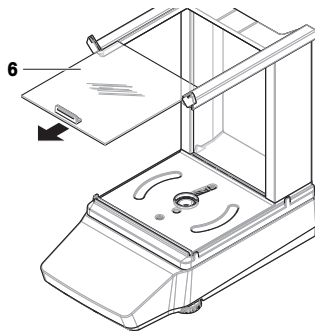
- 5 Pull the top door (6) towards the front to remove it.

- 6 Press the release button (7) and tilt the back panel (8) to remove it.


Note

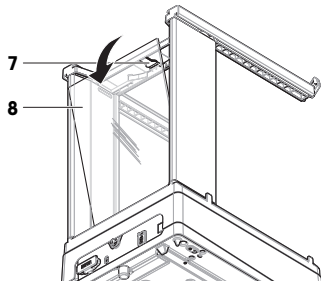Optional, if required: Remove the protective cover for cleaning as described below.


- 7 Remove the front screw (9) and the rear screw (10) with a Phillips screwdriver.

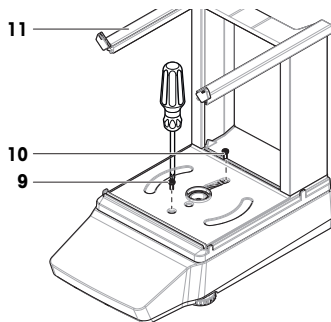- 8 Remove the draft shield (11).

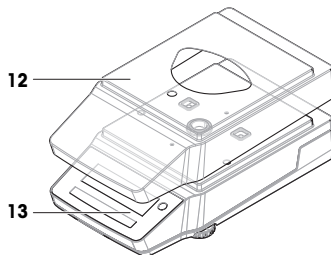- 9 Remove the protective cover (12) from the weighing platform (13).


5.2.1.2 Balances without draft shield


- 1 Remove the draft-protection element (1).
- 2 Remove the weighing pan (2).
- 3 Remove the weighing pan support (3).

Note

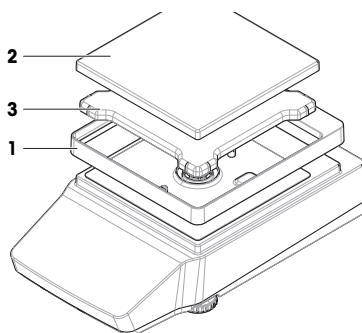Optional, if required: Remove the protective cover for cleaning as described below.

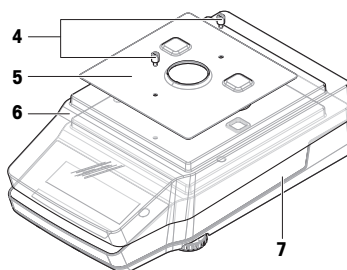
- 4 Remove the screws (4) to remove the EMC plate (5).
- 5 Remove the protective cover (6) from the weighing platform (7).

5.2.1.3 Balances, large

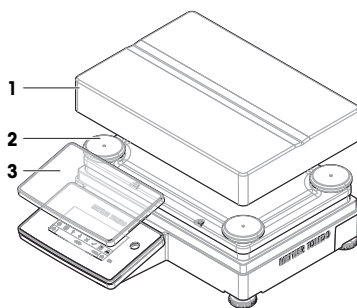- 1 Remove the weighing pan (1).
- 2 Remove the support caps (2).
- 3 Optional, if required: Remove the protective cover (3) for cleaning.


5.2.1.4 Balances, compact


- 1 Remove the weighing pan (1).
- 2 Remove the support caps (2).

Note

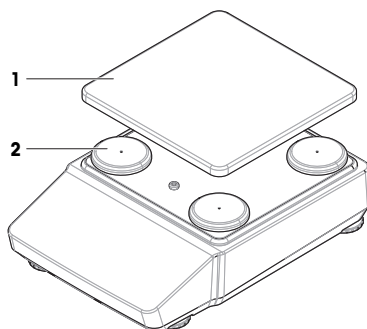Optional, if required: Remove the protective cover for cleaning as described below.

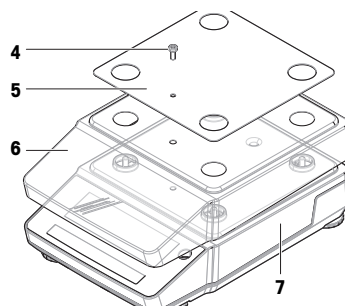- 3 Remove the screw (4) to remove the EMC plate (5).
- 4 Remove the protective cover (6) from the weighing platform (7).

5.2.2 Cleaning the balance

NOTICE

Damage to the instrument due to inappropriate cleaning methods

If liquid enters the housing, it can damage the instrument. The surface of the instrument can be damaged by certain cleaning agents, solvents, or abrasives.

- 1 Do not spray or pour liquid on the instrument.
- 2 Only use the cleaning agents specified in the Reference Manual (RM) of the instrument or the guide "8 Steps to a Clean Balance".
- 3 Only use a lightly moistened, lint-free cloth or a tissue to clean the instrument.
- 4 Wipe off any spills immediately.

For further information on cleaning a balance, consult "8 Steps to a Clean Balance".

► www.mt.com/lab-cleaning-guide

Cleaning around the balance

- Remove any dirt or dust around the balance and avoid further contaminations.

Cleaning the terminal

- Clean the terminal with a damp cloth or a tissue and a mild cleaning agent.

Cleaning the removable parts

- Clean the removed part with a damp cloth or a tissue and a mild cleaning agent or clean in a dishwasher up to 80 °C.

Cleaning the weighing unit

- 1 Disconnect the balance from the AC/DC adapter.
- 2 Use a lint-free cloth moistened with a mild cleaning agent to clean the surface of the balance.
- 3 Remove powder or dust with a disposable tissue first.
- 4 Remove sticky substances with a damp lint-free cloth and a mild solvent, e.g., isopropanol or ethanol 70%.

5.2.3 Putting into operation after cleaning

- 1 Reassemble the balance.
- 2 Check that the draft shield doors (top, sides) open and close normally (if applicable).
- 3 Reconnect the balance to the AC/DC adapter.
- 4 Check the level status, level the balance if necessary.
- 5 Respect the warm-up time specified in the "Technical Data".
- 6 Perform an internal adjustment.
- 7 Perform a routine test according to the internal regulations of your company. METTLER TOLEDO recommends performing a sensitivity test after cleaning the balance.
- 8 Press **→0←** to zero the balance.
 - ➔ The balance is ready for use.

See also

- 🔗 Leveling the balance ▶ Page 19
- 🔗 Technical Data ▶ Page 30
- 🔗 Performing an internal adjustment ▶ Page 21

5.3 Service

Regular servicing by an authorized service technician ensures reliability for years to come. Contact your METTLER TOLEDO representative for details about the available service options.

6 Technical Data

6.1 General data

Power supply

AC/DC adapter:	Input: 100 – 240 V AC ± 10%, 50 – 60 Hz, 0.5 A Output: 12 V DC, 1 A, LPS
Balance power consumption:	12 V DC ± 10%, 0.5 A
Polarity:	

Power supply (optional)

AC/DC adapter:	Input: 100 – 240 V AC ± 10%, 50 – 60 Hz, 0.8 A, 61 – 80 VA Output: 12 V DC, 2.5 A, LPS
Cable for AC/DC adapter:	3-core, with country-specific plug
Balance power consumption:	12 V DC ± 10%, 0.5 A
Polarity:	

Protection and standards

Overvoltage category:	II
-----------------------	----

Degree of pollution:	2
Ingress protection code:	IP43 (precision balances only)

Note

Stated IP is only achieved when the balance is ready for operation. The protective cover must be installed, and the caps must cover the interface connections.

Standards for safety and EMC:	See Declaration of Conformity
Range of application:	Use only indoors in dry locations

Environmental conditions

The limit values apply when the balance is used under the following environmental conditions:

Height above mean sea level:	Up to 5000 m
Ambient temperature:	+10 – +30 °C
Temperature change, max.:	5 °C/h
Relative humidity:	30 – 70%, non-condensing
Acclimatization time:	Recommendation: Up to 4 hours for precision balances, or up to 8 hours for analytical balances. These values apply after placing the balance in the same location where it will be put into operation.

Note

The acclimatization time depends on the readability of the balance, and on the environmental conditions.

Warm-up time:	At least 30 minutes for precision balances, or 60 minutes for analytical balances. These values apply after connecting the balance to the power supply. When switched on from standby, the balance is ready for operation immediately.
---------------	--

The balance can be used under the following environmental conditions. However, the weighing performances of the balance may be outside the limit values:

Ambient temperature:	+5 °C – +40 °C
Relative humidity:	20% to max. 80% at 31 °C, decreasing linearly to 50% at 40 °C, non-condensing

The balance can be disconnected and stored in its packaging under the following conditions:

Ambient temperature:	-25 – +70 °C
Relative humidity:	10 – 90%, non-condensing

7 Disposal

In conformance with the European Directive 2012/19/EU on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device. Should this device be passed on to other parties, the content of this regulation must also be related.

8 Compliance Information

National approval documents, e.g., the FCC Supplier Declaration of Conformity, are available online and/or included in the packaging.

► www.mt.com/ComplianceSearch

For further information, consult the Reference Manual (RM).

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► www.mt.com/contact

1.3 Akronymmer og forkortelser

Originalt udtryk	Oversat udtryk	Forklaring
AC		Alternating Current
ASTM		American Society for Testing and Materials
DC		Direct Current
EMC		Electromagnetic Compatibility
FCC		Federal Communications Commission
GWP		Good Weighing Practice
HID		Human Interaction Device
ID		Identification
IP		Ingress Protection
LAN		Local Area Network
LED		Light-Emitting Diode
LPS		Limited Power Source

MAC	Media Access Control
MT-SICS	METTLER TOLEDO Standard Interface Command Set
NA	Not Applicable
OIML	Organisation Internationale de Métrologie Légale (International Organization of Legal Metrology)
RAM	Random Access Memory
RM	Reference Manual
SOP	Standard Operating Procedure
UM	User Manual (Brugervejledning)
USB	Universal Serial Bus
USP	United States Pharmacopeia

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© 12/2023 METTLER TOLEDO. All rights reserved.